



WRO 300 INLET HOSE MODIFICATION

The WRO 300/300 H unit is supplied with an installation kit, W3T578238 or W3T579082 depending on model. This kit includes a quick disconnect hose barb fitting. The cart contains an inlet supply hose, PN W3T577916. This hose needs to be modified and quick connect fitting attached as instructed below. This Tech Note applies to M1, M2 Ergo and Ergo Docking Station

1. To connect the outlet of the pretreatment filter to the inlet of the WRO 300 requires modification of the hose W3T577916.

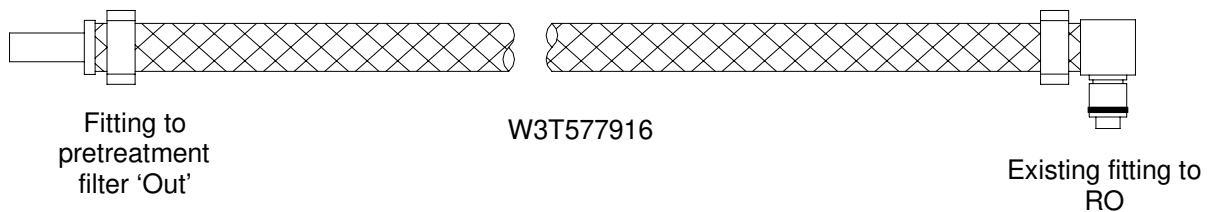

2. Measure approximately 6" from the pretreatment filter fitting end of hose. Cut hose. Keep the 6" length of hose with the pretreatment fitting end. Discard the remaining length of hose with the existing QD (quick disconnect).

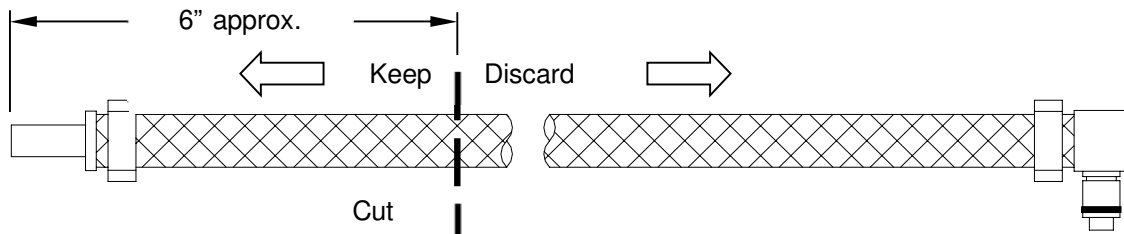

3. Install straight male QD from WRO 300 install kit using hose clamp from same kit.

