

### **WRO 300 / WRO 300 H WATER SYSTEMS TIME CHANNEL PROGRAM #14**

#### **Purpose**

The purpose of this Tech Note is to inform users of an issue with Time Channel Program #14.

Time Channel Program #14 (GXP parameter S215) in the WRO 300 and WRO 300 H water systems cannot be used due to a software code typo that prevents it from being saved. The time channel can be setup for the various operating modes and times, however, when saving TO time channel #14 it will indicate an error has occurred.

The remaining 20-time channels (1-13 and 15-21; S159-S165, S209-S214 and S216-S222) remain fully functional. This issue occurs when setting time channels through the user interface or when using the GXP software. There are no other impacts to the software for the WRO 300 and WRO 300 H Water Purification Systems.